


Upgrading Stereotaxic Instruments -2

A great variety of upgrades are available for our stereotaxic instruments.
In this issue, we introduce new optional accessories.

◆◇ For Gas Anesthesia ◆◇

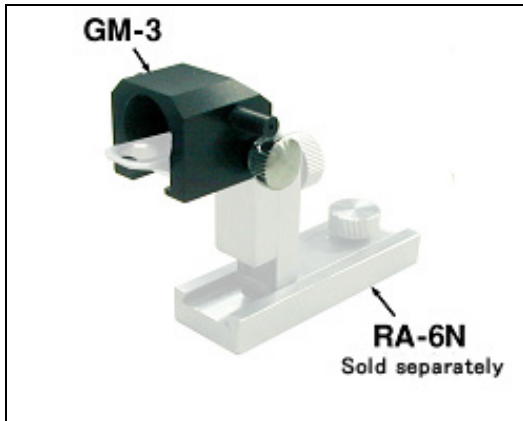


When an animal should go under anesthetic state in an experiment, you can use the **GM-3** (for rats) or **GM-4** (for mice) to connect with 4mm I.D. anesthetic tube to permit gas anesthesia. As the anesthetic mask holds the nose by itself, three-point fixation is maintained to hold the head solidly.

*MRI-compatible anesthetic mask is also available as special order.

◆◇ For Cannula Placement ◆◇

The **SHC-1** is designed to be mounted on micromanipulator on stereotaxic instrument to allow quick and accurate placement of cannula without annoying hand vibration. It is also designed to allow easy holding and release of a cannula with single knob, and accept almost all 3.5mm diameter cannulas available on the market. We can also produce cannula holders for other specifications as special order. If you have any questions, please feel free to consult us.

*We are happy to advise on compatibility. When you inquire, please tell us cannula model.

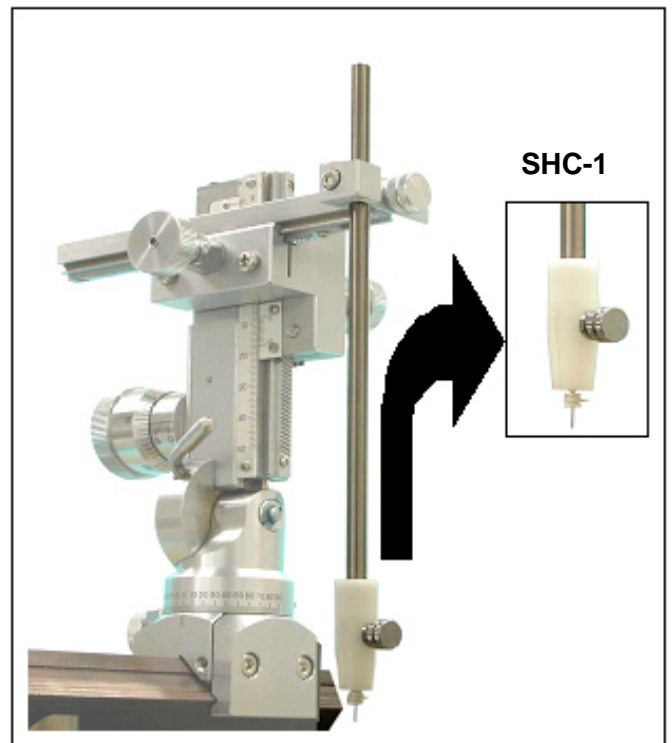

☆☆ What is cannula? ☆☆

It is a tube which is used to inject drug or electrode into brain, or collect body fluid in animal testing.